Vigtig produktinformation

KORREKTION nr. 1218959-10-18-2012-006-C

Rettelse til medicinsk enhed

18. oktober 2012

Bemærk: Kunde

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Waters ACQUITY® TQD, ACQUITY® SQD, LCT Premier™ XE, Quattro Premier™, Quattro Premier™ XE, Quattro™ mikro, Xevo® G2 Tof, Xevo® TQ, Xevo® TQD & ZQ 2000 massespektrometre, alle serienumre.

Beskrivelse af problemet:

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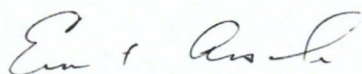
Venlig hilsen

Ernest Ansah
Quality Assurance Manager, Clinical Quality
Waters Corporation
34 Maple Street
Milford, MA 01757
Telefon: +1 508-482-2811
Fax: +1 508-482-2339

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Description of the problem:

Waters Corporation is initiating a voluntary medical device correction for its mass spectrometers referenced above. Our records indicate that you own or lease one or more of these mass spectrometers.

This correction is necessary because there is a possibility that the ion block source heater may fail. In some instances, the failure could cause a short term loss of or variation in sensitivity that may subsequently lead to inaccurate data. The loss of or variation in sensitivity is most pronounced in assays with shorter dwell times (i.e. less than 100 ms) and may not always be detected by quality control and internal standard value monitoring.

Advise on action to be taken by the user:

To ensure the continued safety and reliability of our products Waters recommends that you replace the ion block source heater immediately. The necessary part(s) are included with this Field Safety Notice and Waters estimates the replacement should take approximately 30 minutes. The ion block source heater is customer replaceable as detailed in the maintenance procedure of the operators guide. Instructions can be also be found in several languages at the following address: www.waters.com/sourceheater. Alternatively, a local field service engineer can replace the ion block source heater at no cost. If you prefer this alternative, please contact your local field service support office¹ to schedule this service.

Please complete the attached Reply Verification Tracking Form and return it to Waters via e-mail at waters_quality@waters.com, or via fax at the number provided below. This form provides documented evidence that this notification and the replacement part were received at your site.

¹ To contact your local field service engineer outside of the United States, log onto the Waters Website at <http://www.waters.com/waters/home.htm>; choose your country at the top of the page, click on About Waters/Contact Waters/Local Country Contacts.

Transmission of this Field Safety Notice:

This notice needs to be passed on to all those who need to be aware within your organisation or to any organisation where the potentially affected instruments have been transferred.

Contact reference person:

The undersigned confirms that this notice has been provided to the appropriate Regulatory Agency.

We regret any inconvenience that this action may cause. Please do not hesitate to contact me if you have any questions.

Thank you,

Ernest Ansah
Quality Assurance Manager, Clinical Quality
Waters Corporation
34 Maple Street
Milford, MA 01757
Telephone: +1 508-482-2811
Fax: +1 508-482-2339

**Medical Device
Correction**

REPLY VERIFICATION TRACKING FORM

Re: Waters ACQUITY® TQD, ACQUITY® SQD, LCT Premier™ XE, Quattro Premier™, Quattro Premier™ XE, Quattro™ micro, Xevo® G2 Tof, Xevo® TQ, Xevo® TQD & ZQ 2000 Mass Spectrometers, all serial numbers.

INSTRUCTIONS FOR COMPLETING THIS REPLY VERIFICATION TRACKING FORM:

This completed form is required for tracking purposes. **Please check the applicable boxes, provide the required information and return this form** via Fax or email to:

- (f) Fax Number: (508) 482-2339 or
- (e) E-mail address: waters_quality@waters.com

SECTION TO BE COMPLETED BY CUSTOMER

- We received the notification and the replacement ion block source heater
- The ion block source heater was replaced

Company Name _____

Company Address _____

Contact name: _____ Other: _____

Contact Title: _____

Department: _____ Telephone Number: _____

Signature (required): _____ Date: _____

Email address: _____

Reply Verification Tracking Form Serial Number: _____

Field Action Identification Number 1218959-10-18-2012-006-C

120F12